

Laser Speckle Reducer LSR-3000 Series



Speckle noise from a laser-based system is reduced by dynamically diffusing the laser beam. A diffuser is bonded to a thin elastic membrane, which includes four independent electro-active polymer electrodes that induce a circular oscillation of the diffuser in x- and y-direction. The LSR-3000 Series integrates fully certified drive electronics powered through a single micro-USB connector. Two sizes are available: the LSR-3005 and the LSR-3010 exhibiting a clear aperture of 5 mm and 10 mm, respectively.

Mechanical specifications

	LSR-3005	LSR-3010	
Clear aperture	5	10	mm
Outer diameter	41	48	mm
Thickness	8.8	8.8	mm
Weight	24.43	32.84	g

Electrical specifications¹

Power supply (micro-USB interface)	5	5	VDC
Power consumption	310	310	mW

Optical specifications

Diffusion angle (FWHM)	6°/12°/17°/24°	6°/12°/17°/24°	
Transmission spectrum ²	see figures 2&3	see figures 2&3	
Damage threshold	>300	>300	W/cm ²
Oscillation frequency	~300	~180	Hz
Oscillation amplitude (peak to peak)	~300	~400	µm

Environmental specifications

Storage temperature	[-40,+85]	[-40,+85]	°C
Operating temperature	[-30,+85]	[-30,+85]	°C
Lifetime in operation ²	2000	2000	h

Figure 1 (a-c) show the effect of the LSR-3005 on a laser spot ($\lambda = 650$ nm, $P = 5$ mW).

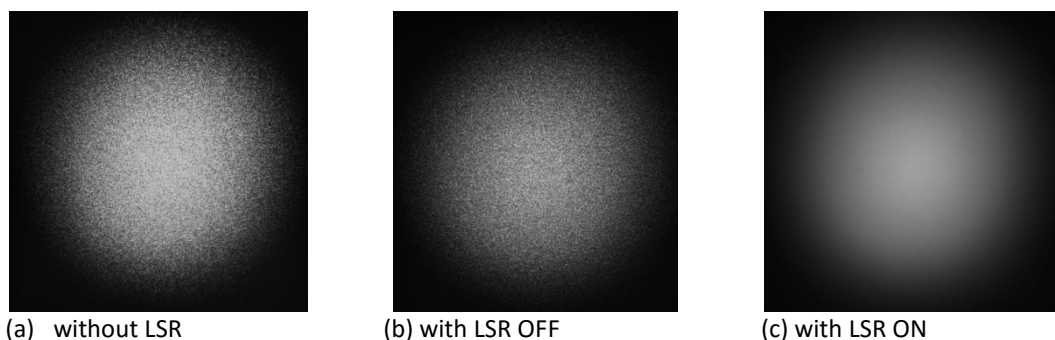


Figure 1: CCD images of a laser spot without and with LSR-3005 on.

¹ 100-230 VAC to 5 VDC micro-USB power supply provided by Optotune

² This product is ideal for applications with a low duty cycle. For applications with long lifetime Optotune's reluctance force LSR products are recommended

The following two figures show the transmission spectrum of the LSR-3000 series with polymer diffusers and coated BK7 cover glasses. The LSR-3000-xxD series include both a static and an oscillating diffuser for increased speckle reduction efficiency. Only an oscillating diffuser and no static diffuser are present in the LSR-3000-xxS series.

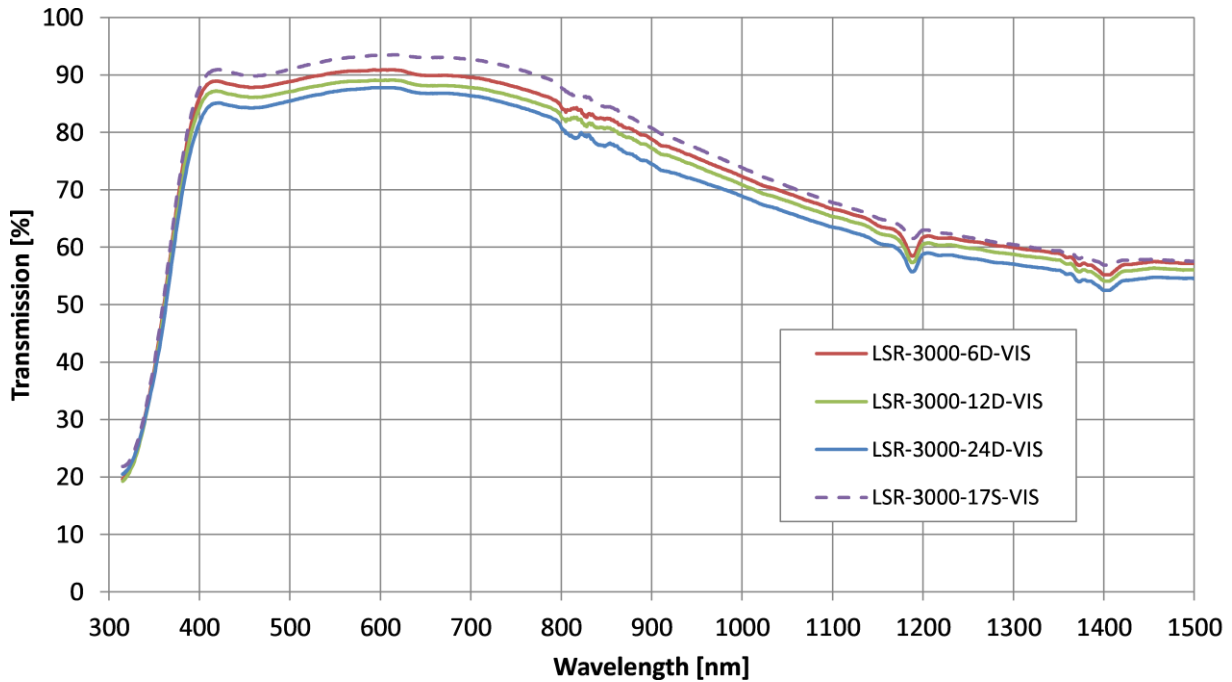


Figure 2: Transmission spectrum of the LSR-3000 series with VIS-coated cover glasses (0° angle of incidence)

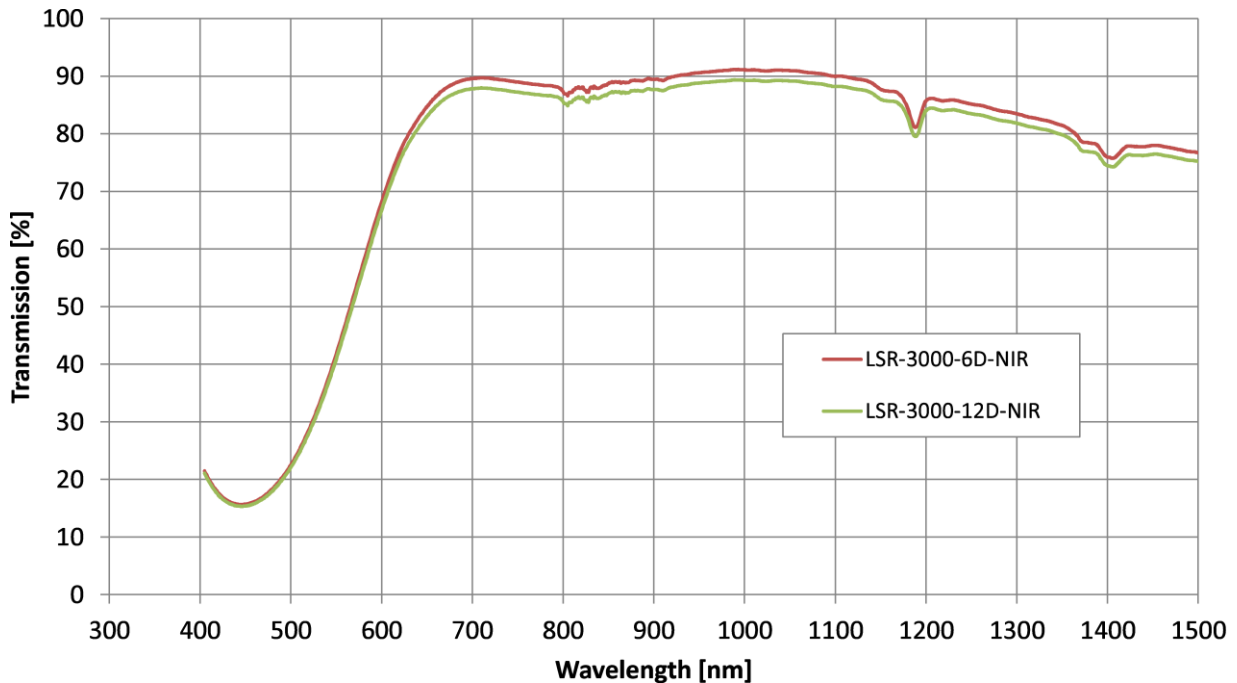


Figure 3: Transmission spectrum of the LSR-3000 series with NIR-coated cover glasses (0° angle of incidence)

Figure 4 and Figure 5 show the dimensions of the LSR-3005 and the LSR-3010, respectively. The housing exhibits a M4 thread, in order to facilitate the integration of the LSR on an optical table. The power supply interface is made through a micro-USB connector. The LSR-3000 series is also available without housing (see datasheet of LSR-OEM).

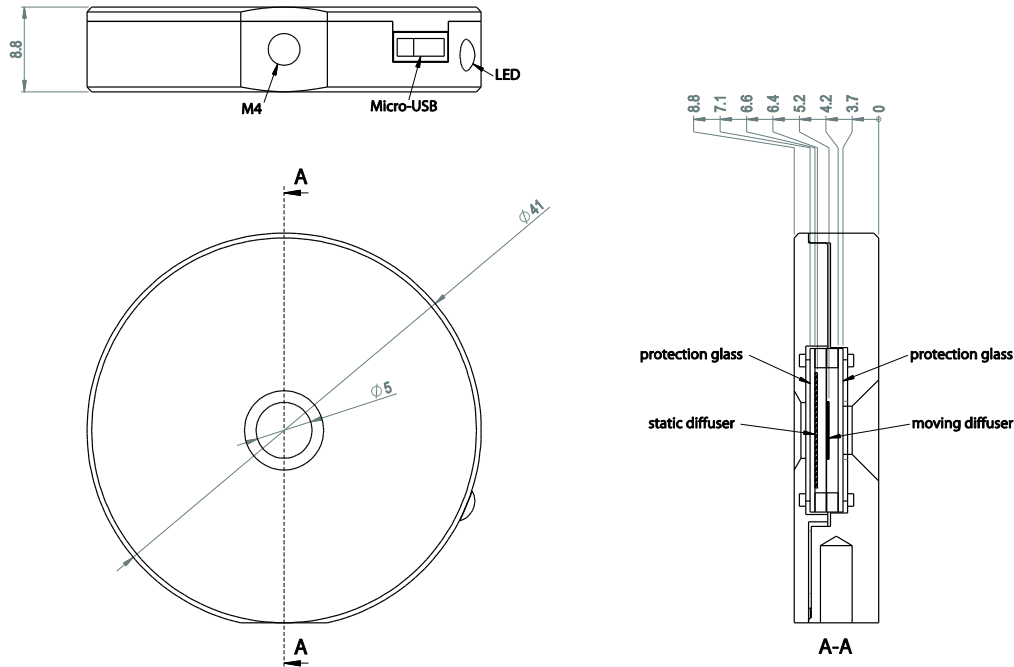


Figure 4: Mechanical drawing of the LSR-3005 (unit: mm)

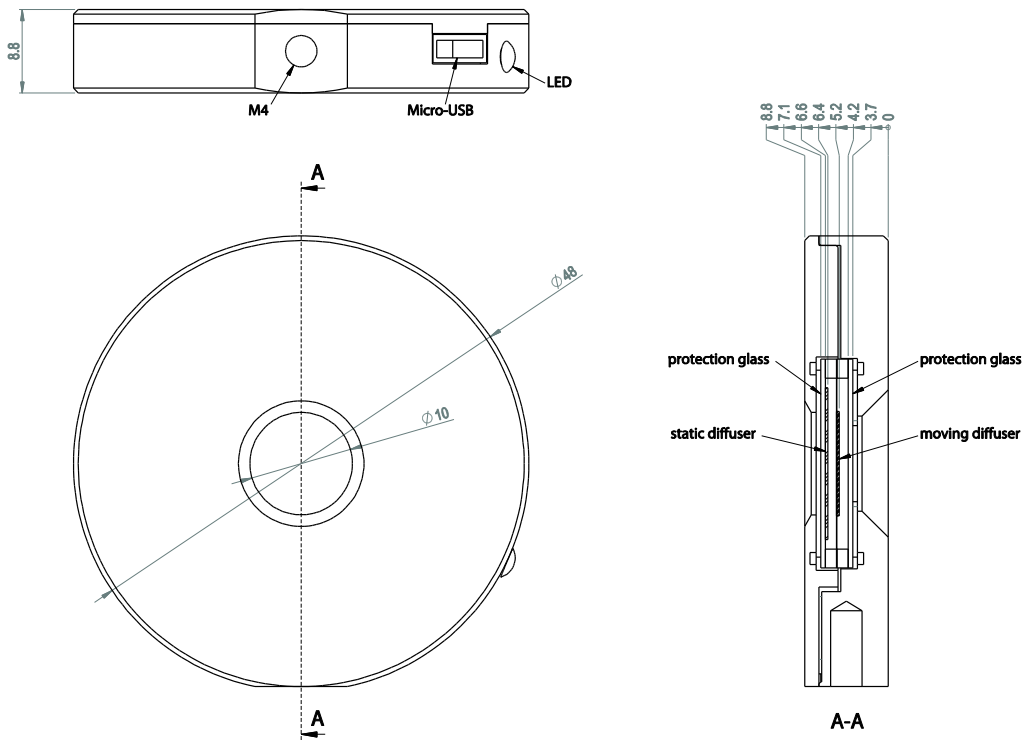


Figure 5: Mechanical drawing of the LSR-3010 (unit: mm)

Lifetime

The LSR has passed the following accelerated aging tests. The tests were performed on 5-10 samples each and are considered as passed if the oscillation amplitude of all samples did not change more than 5%.

Test	LSR-3005
A1 – High temperature (operating): <ul style="list-style-type: none"> Test temperature: 85°C ±5°C Test duration: 48 hours Reference: JESD22-A108-B, JESD47-A and JESD74 	Passed
A2 – Low temperature (operating): <ul style="list-style-type: none"> Test temperature: -30°C ±5°C Test duration: 48 hours Reference: JESD22-A108-B, JESD47-A 	Passed
A3.1 – Temperature cycling (non-operating): <ul style="list-style-type: none"> Low temperature: -40°C, high temperature: +85°C Transition time: <3 minutes Test duration: 30 min for each under low and high temperature Number of cycles: 100 cycles Recovery time: 2 hours References: IEC 60068-2-1, IEC 60068-2-2 and IEC 60068-2-14 	Passed
A3.2 – Damp heat cycling (non-operating): <ul style="list-style-type: none"> Low temperature: 25°C, high temperature: +55°C Test humidity: 90-100% relative humidity Test duration: 12 hours dwell Temperature change rate: Less than 1°C per minute Number of cycles: 18 cycles Recovery time: No recovery time References: IEC 60068-2-30 Db Variant 1 	Passed
A4 – High temperature (non-operating): <ul style="list-style-type: none"> Test temperature: Static 85°C ±2°C Test humidity: Less than 20 g/m³ of water vapor (corresponding to approximately 50% relative humidity at 35°C) Test duration: 168 hours Temperature change rate: Less than 1°C per minute Reference: IEC 60068-2-2 Bb 	Passed
Shock test: <ul style="list-style-type: none"> 400 g for 1 ms duration 5 pulses in each direction (30 pulses in total) 	Passed
Operating lifetime test: <ul style="list-style-type: none"> Test temperature: 22 ±2°C Test duration: 2000 hours Driving frequency: 300 Hz 	Passed ³

Table 1: Environmental and accelerated aging tests performed on the LSR Series

³ The average amplitude of 300um dropped to 200um after 2000 hours of operation.

Ordering information for standard products

The LSR-3000 can be delivered with two sizes of clear aperture, different diffusing angles (circular) and two types of cover glasses. When ordering, please refer to these specifications as follows:

LSR-30CA-θN-AR

CA = Clear aperture (05 or 10 mm)

θ = Total diffusion angle (6°, 12°, 17°, 24° available as standard)

N = Diffuser configuration, whereas D = double (one oscillating, one static), S = single (one oscillating only)

AR = Antireflection coating (VIS or NIR)

Example: LSR-3005-12D-VIS refers to a speckle reducer of 5-mm aperture with a total diffusion angle of 12°, with both an oscillating and a static diffuser and VIS-coated cover glasses.

Standard models with different diffuser combinations

The standard models combine an oscillating diffuser with a static diffuser to reduce the correlation length of the random patterns that are generated. This provides more effective speckle reduction yet minimizing the increase in beam divergence. Three standard diffuser configurations are available: LSR-3000-xxD series with 6° (FWHM) for highest optical efficiency and 24° for best speckle reduction. For optical systems where the spot on the LSR is imaged e.g. by a 4f-system, no static diffuser is allowed. In that case, the LSR-3000-xxS series with a oscillating diffuser only is recommended.

Standard products	Total diffusion angle	Diffuser configuration	Cover glass coating
LSR-3010-6D-VIS	6°	4.2° oscillating, 4.2° static	400 – 700 nm
LSR-3010-12D-VIS	12°	8.5° oscillating, 8.5° static	400 – 700 nm
LSR-3005-6D-VIS	6°	4.2° oscillating, 4.2° static	400 – 700 nm
LSR-3005-12D-VIS	12°	8.5° oscillating, 8.5° static	400 – 700 nm
LSR-3005-24D-VIS	24°	17° oscillating, 17° static	400 – 700 nm
LSR-3005-17S-VIS	17°	17° oscillating, no static	400 – 700 nm
LSR-3005-6D-NIR	6°	4.2° oscillating, 4.2° static	700 – 1100 nm
LSR-3005-12D-NIR	12°	8.5° oscillating, 8.5° static	700 – 1100 nm

Table 2: Overview of standard LSR-3000 models

Custom products

Optotune offers customized versions of the LSR. This datasheet only contains variations of the LSR-3000 Series products. For LSRs with substantially smaller or larger apertures, please contact sales@optotune.com.

Diffusers

In principle, any type of diffuser can be used as long as size and weight are similar to the standard diffuser. Furthermore, it is possible to use specialty coated glass diffusers if they are light enough. If necessary, it is also possible to remove the membrane from the optical path.

The following table summarizes these options:

Diffuser type	Optotune polymer	Glass diffusers (typical)
Thickness	330 µm	As thin as possible
Density	1.03 g/cm ³	2.2 g/cm ³
Weight		
5 mm aperture ⁴	9 mg	up to 20 mg
10 mm aperture	30 mg	up to 70 mg
Resonance frequency		
5 mm aperture	300 Hz	$\omega \propto \sqrt{\frac{1}{m}}$
10 mm aperture	180 Hz	
Transmission range	240 – 2500 nm	170 – 12000 nm (depending on material)
Damage threshold	> 300 W/cm ²	1 kW/cm ² (depending on material)

Please note that diffusers cannot be exchanged after the LSR has been produced. Also, when using customized diffusers it might be necessary to select an appropriate cover glass to match the transmission range.

Cover glasses

Cover glasses can be customized with respect to material and AR-coating. The standard cover glasses are VIS or NIR coated BK7. If the LSR is to be integrated into a clean environment, then the cover glasses can be omitted.

Frequency

The oscillation frequency is set to the measured resonant frequency of the LSR during production. With the current standard electronics it is not possible to change that frequency after production.

Safety and compliance

The product fulfills the RoHS and REACH compliance standards. The customer is solely responsible to comply with all relevant safety regulations for integration and operation.

For more information on optical, mechanical and electrical parameters, please contact sales@optotune.com.

⁴ Note: The 5mm and 10mm aperture LSR is usually equipped with 6mm and 11mm diameter diffuser, respectively.